EK-*i* / EW-*i*

Quick Start Guide
EK-*i* / EW-*i* Series
Compact Balances

English (EN)

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The logo for A&D Company, Ltd. features the letters 'A', 'D', and '&' in a bold, blue, sans-serif font. The 'A' and 'D' are significantly larger than the ampersand, and they are all connected together.

A&D Company, Ltd.

1WMPD4004399

Quick Start Guide – EK- <i>i</i> / EW- <i>i</i> Series	EN
Compact Balance	3
Guide de démarrage rapide – Séries EK- <i>i</i> / EW- <i>i</i>	FR
Balance compactes	12
Schnellstartanleitung – Serie EK- <i>i</i> / EW- <i>i</i>	DE
Kompakte-Waage	21
Guía de inicio rápido – Series EK- <i>i</i> / EW- <i>i</i>	ES
Básculas compactas	30
Guida introduttiva – Serie EK- <i>i</i> / EW- <i>i</i>	IT
Bilancia compatte	39
快速入門指南- EK- <i>i</i> / EW- <i>i</i> 系列	CT
緊湊型天平	48

1. Introduction

This quick start guide describes the basic functions and operations of the balance.

The functions and operating procedures of the balance (EK-*i* / EW-*i* series) are explained in details in the following instruction manual.

[EK-*i* / EW-*i* Series Instruction Manual]

This is an instruction manual for understanding and fully utilizing the precision electronic balance EK-*i* / EW-*i* series. This manual can be downloaded from our website <https://www.aandd.jp>.


2. Unpacking

When unpacking, check whether all of the following items are included:

Weighing pan

Pan support


Main unit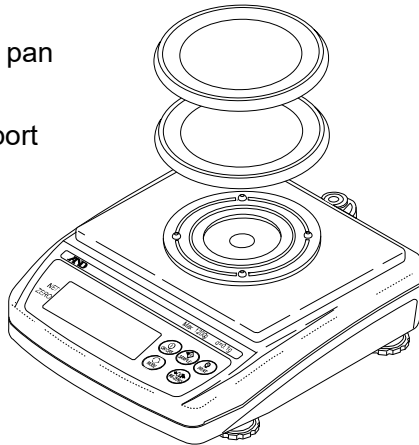

EK-120*i* / EK-200*i* / EK-300*i*
 EK-410*i* / EK-610*i*
 EW-150*i*


Weighing pan

Main unit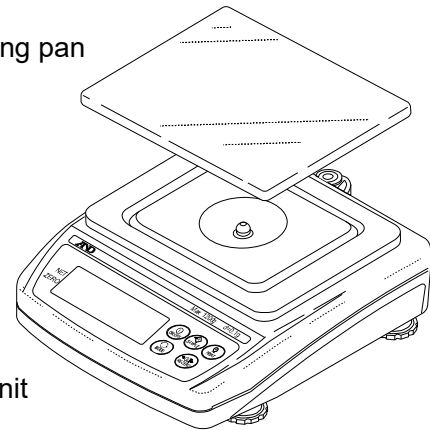

EK-600*i* / EK-1200*i* / EK-2000*i*
 EK-3000*i* / EK-6000*i* / EK-12K*i*
 EK-4100*i* / EK-6100*i*
 EW-1500*i* / EW-12K*i*


AC adapter

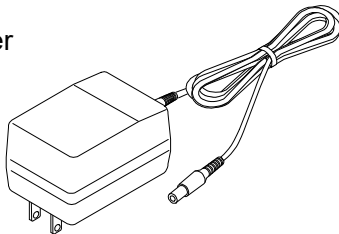
Note:

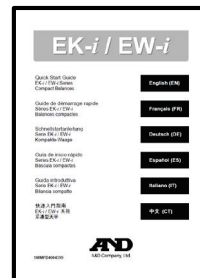Verify that the AC adapter type is correct for your local voltage and receptacle type.

Quick Start Guide

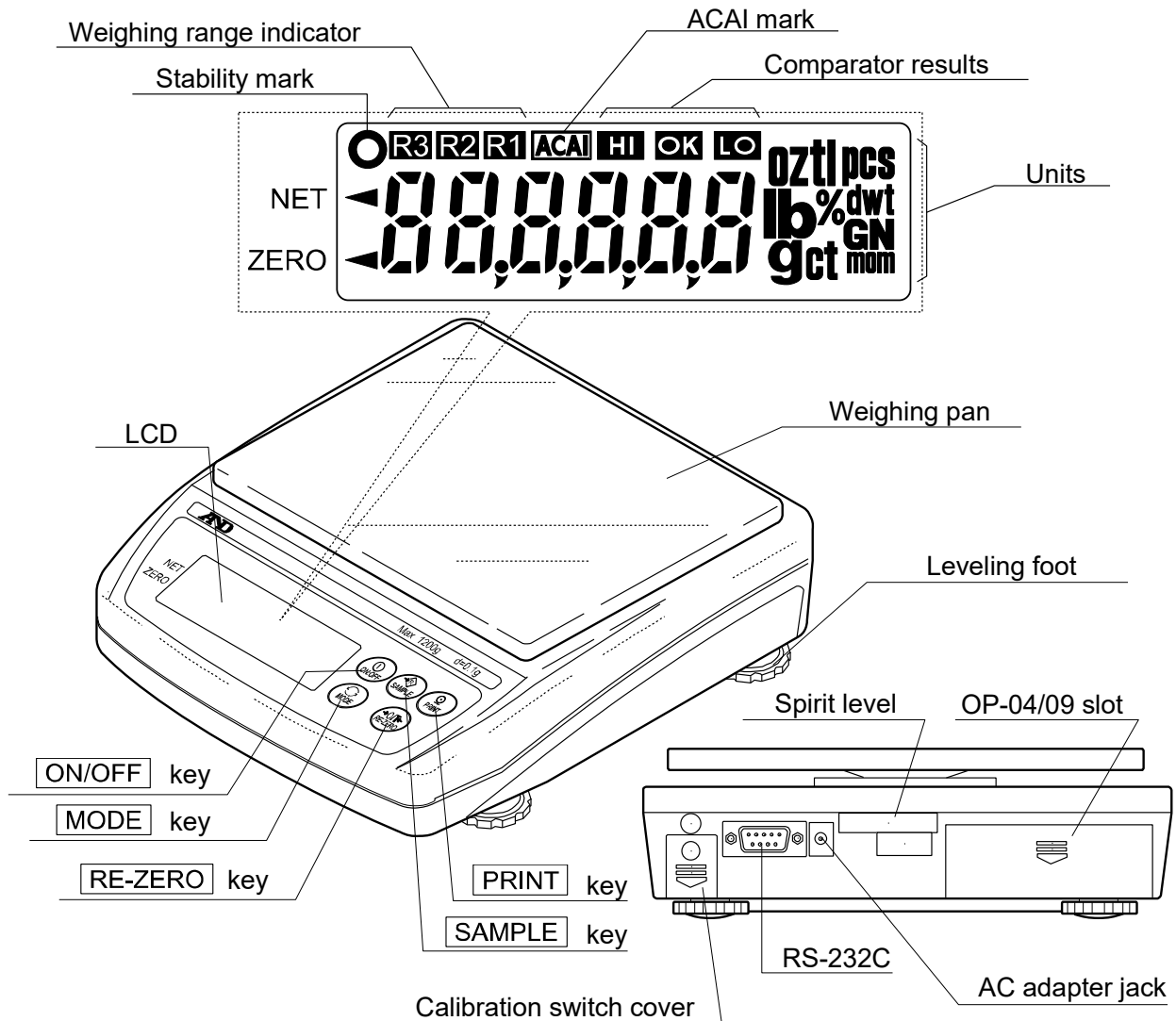3. Part names and functions

 <p>Turns the power ON or OFF.</p>	 <p>Outputs the weight data to a printer or computer, or enters a unit weight, 100% weight or other setting values to the balance.</p>
 <p>Hold down to enter the function setting mode.</p> <p>Weighing mode (EW-<i>i</i> only): Changes the weighing range (when rnG 0 is selected).</p> <p>pcs mode: Enters the sample unit weight storing mode.</p> <p>% mode: Enters the 100% weight storing mode.</p>	 <p>Switches the weighing unit (the weighing mode).</p>  <p>Clears the display to zero.</p>

4. Setting up

4-1 Setting up your balance

1. Place the weighing pan on the main unit as shown on the previous page.
2. Adjust the level of the balance using the leveling feet.
Use the spirit level to confirm. The bubble should be in the center of the circle.
3. Calibrate your balance before use. (See “6. Calibration”)

Balance installation

To measure correctly, to keep the balance in good condition, and to prevent hazards, observe the following:

- Do not install the balance in locations that are subject to excessive dust, breezes, vibration, large temperature fluctuations, condensation, or that may have magnetic fields.
- Do not install the balance on a surface that is soft or that may cause the balance level to shift.
- Do not install the balance in direct sunshine.
- Do not install the balance near heaters or air conditioners.
- Do not use an unstable AC power source.
- Do not install the balance in a place where combustible or corrosive gases may exist.
- Allow the balance to reach equilibrium with the ambient temperature before use.
- Switch the power ON at least half an hour before use so that the balance can warm up.
- When the balance is installed for the first time, or the balance has been moved, carry out calibration as described in “6. Calibration”.

4-2 Power source

For the power source, the AC adapter or the rechargeable battery pack (EKW-09*i*: Optional item) is available.

When using the AC adapter

Use a stable power source. To use the AC adapter, insert the AC adapter plug into the AC adapter jack on the EK-*i* / EW-*i*.


When using the rechargeable battery pack (EKW-09*i*)

Refer to the [EK-*i* / EW-*i* Series Instruction Manual] for more information.

5. Operation

5-1 Turning the power ON and OFF

1. Press the key to turn the power ON.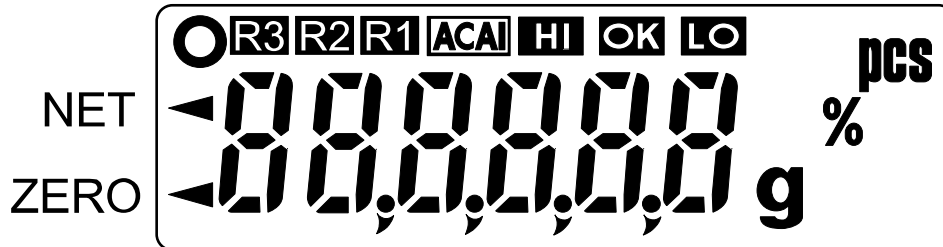

All of the indicators are displayed as shown above.

(About units: Only the units available will be displayed.)

The display turns off except for a weighing unit and the decimal point.

The balance waits for the weight data to become stable, and zero will be displayed with the ZERO mark (power-on zero).

The range for power-on zero is within $\pm 10\%$ of the weighing capacity around the calibrated zero point.

If the power is switched ON while there is a load beyond this range, the balance will be tared to zero and the NET mark and the ZERO mark turn on.

2. Pressing the key again, and the power will be switched OFF.


6. Calibration

This function adjusts the balance for accurate weighing.

Perform calibration in the following cases.

- When the balance is first used.
- When the balance has been moved.
- When the ambient environment has changed.
- For regular calibration.

Press and lower the calibration switch cover.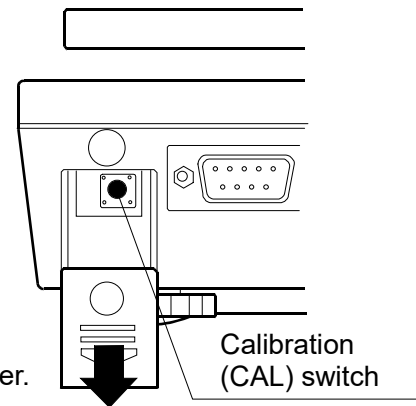

6-1 Calibration using a weight

- Prepare a calibration weight (optional) before start.
1. Warm up the balance for at least half an hour with nothing on the pan.

2. Press and hold the calibration (CAL) switch until **[CAL]** appears, and release the switch.

3. The balance displays **[CAL 0]**.

To change the calibration weight value, proceed to step 4.

To use the calibration weight value in the balance memory, proceed to step 5.

4. Press the **[SAMPLE]** key. The display shows the calibration weight value in “gram” that is stored in the balance. Use the following keys to change the value.

[SAMPLE] key To select the digit blinking to change.

[RE-ZERO] key To set the value of the digit selected.

[PRINT] key To store the value and return to step 3.

[MODE] key To cancel the value and return to step 3.


- Use a calibration weight of more than 2/3 of the capacity (of the highest range for EW- *i* series).

 Press and hold the CAL switch.

[CAL]


Release the CAL switch.

[CAL 0]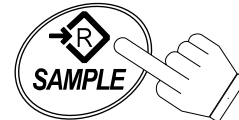

[0 12000]

Set the weight using the relevant keys.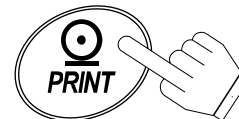


[CAL 0]


5. At step 3, pressing the **PRINT** key weighs the zero-point value. Do not touch the pan during weighing.

When the zero calibration is completed, the display shows the calibration weight value.

- To perform the zero calibration only and finish the procedure, press the **MODE** key.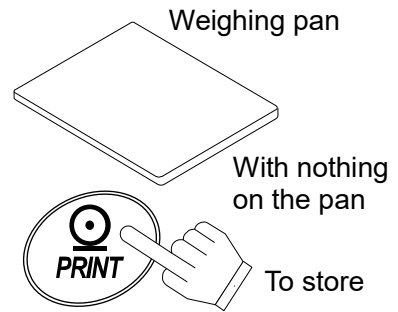

6. Place the calibration weight with the same value as displayed on the pan. Press the **PRINT** key to weigh it. Do not touch the pan during weighing.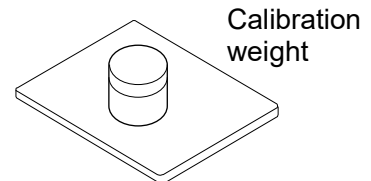

7. **End** appears.

Remove the weight from the pan, and press the CAL switch or **MODE** key to return to the weighing mode.










7. Maintenance

7-1 Notes on maintenance

- Do not disassemble the balance. Contact your local A&D dealer if your balance needs service or repair.
- Use the original package for transportation.
- Do not use organic solvents to clean the balance. Use a lint free cloth dampened with a mild detergent.

7-2 Error codes

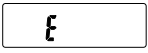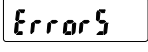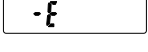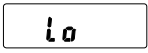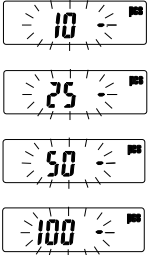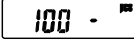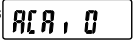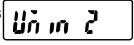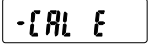Display	Description
	<p>Overload error</p> <p>Warning to indicate that an object beyond the balance capacity has been placed on the pan. Remove the object from the pan.</p>
	<p>Overload error</p> <p>Warning to indicate that the balance detects an unexpected load. Remove all objects from the pan or check that the weighing pan is correctly installed.</p>
	<p>Range over notice</p> <p>This will be shown if the weight sensor receives strong force upward. Check if there is anything sandwiched around the weighing pan. There is a possibility that the weight sensor itself may have a failure.</p>
	<p>Unit / 100% weight error</p> <p>The unit / 100% weight is too light in the counting / % mode.</p>
	<p>Sample quantity notice</p> <p>When sample weight is light and the counting error could become large, the balance will prompt you to use a larger number of samples. Place the displayed number of samples on the pan and press the PRINT key to store the unit weight.</p> <p>Note:</p> <p>Pressing the PRINT key without adding samples may reduce counting accuracy. Starting from the 100 samples,  may be displayed when the sample weight is light. This is for your notice and press the PRINT key without adding any samples.</p> <p>This notice is not shown when " (ACAI disabled) or " is set.</p>
	<p>CAL errors</p> <p>Warning to indicate that calibration has been canceled because the calibration weight is too light.</p> <p>Check the weighing pan and the calibration weight.</p> <p>To return to the weighing mode, press the MODE key.</p>

Lb0	<p>Low battery</p> <p>Warning to indicate that the battery pack (OP-09) is discharged. Immediately stop using it, and recharge or use the AC adapter.</p>
Hb1	<p>AC adapter error</p> <p>Warning to indicate that the output voltage of the AC adapter is too high. Check if the AC adapter is correct.</p>
Lb1	<p>AC adapter error</p> <p>Warning to indicate that the output voltage of the AC adapter is too low. Check if the AC adapter is correct.</p>
Error1	<p>Stability error</p> <p>Warning to indicate that the weight value is not stable and the balance cannot display it. Prevent vibration and drafts. Press the MODE key to return to the weighing mode.</p>
Error#	<p>Internal errors (# = 2, 3, 4 or 6)</p> <p>The balance detects error state in the internal processing. With nothing on the weighing pan, turn the power OFF and ON again. If the error persists, request service.</p>

If you cannot cancel an error or other errors have occurred, request service from the store where you purchased the balance or from your local A&D dealer.

8. Specifications

8-1 EK-*i* series

Model	EK-120 <i>i</i>	EK-200 <i>i</i>	EK-300 <i>i</i>	EK-600 <i>i</i>
Weight capacity	120g	200g	300g	600g
Readability	0.01g	0.01g	0.01g	0.1g
Repeatability (Standard deviation)	0.01g	0.01g	0.01g	0.1g
Linearity	±0.01g	±0.01g	±0.02g	±0.1g

Model	EK-1200 <i>i</i>	EK-2000 <i>i</i>	EK-3000 <i>i</i>	EK-6000 <i>i</i>	EK-12K <i>i</i>
Weight capacity	1200g	2000g	3000g	6000g	12kg
Readability	0.1g	0.1g	0.1g	1g	1g
Repeatability (Standard deviation)	0.1g	0.1g	0.1g	1g	1g
Linearity	±0.1g	±0.1g	±0.2g	±1g	±1g

Model	EK-410 <i>i</i>	EK-610 <i>i</i>	EK-4100 <i>i</i>	EK-6100 <i>i</i>
Weight capacity	400g	600g	4000g	6000g
Readability	0.01g	0.01g	0.1g	0.1g
Repeatability (Standard deviation)	0.01g	0.01g	0.1g	0.1g
Linearity	±0.02g	±0.02g	±0.2g	±0.2g

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A&D Company, Limited
1756 Automation Parkway, San Jose, California 95131, U.S.A.
A&D AUSTRALASIA PTY LTD
A&D KOREA Limited