FG-CWP Series

Quick Start Guide
FG-CWP Series
For Water proof platform scale

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A&D Company, Ltd.

© 2022 A&D Company, Limited. All rights reserved.

No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company, Limited.

The contents of this publication and the specifications of the instrument covered by this publication are subject to change for improvement without notice.

Quick Start Guide – FG-CWP Series	EN
Water proof digital platform scale.....	4
Guide de démarrage rapide – Séries FG-CWP	FR
Balance numérique à plateforme numérique étanche.....	11
Schnellstartanleitung – Serie FG-CWP	DE
Wasserdichten digitalen Plattformwaage.....	19
Guía de inicio rápido – Series FG-CWP	ES
Báscula de plataforma digital impermeable.....	27
Guida introduttiva – Serie FG-CWP	IT
Bilancia digitale impermeabile a piattaforma.....	35
快速入門指南- FG-CWP 系列	CT
防水數位平台秤.....	43

1. Introduction

This quick start guide describes the basic functions and operations of the balance.

The functions and operating procedures of the scale (FG-CWP series) are explained in details in the following instruction manual.

[FG-CWP Series Instruction Manual]

(<https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html>)

This is an instruction manual for understanding and fully utilizing the water proof digital platform scale FG-CWP series. This manual can be downloaded from our website <https://www.aandd.jp>.

2. Cautions

2.1. Precautions for Installing the Scale

DANGER

- Ground the scale, so that the user will not be subjected to an electric shock.
- To prevent electric shock, do not handle the USB cable with wet hands.
- The USB cable is not waterproof. Install the scale in an area where it will not get wet.
- Do not install the scale where flammable or corrosive gas is present.
- Do not put excessive force on cables.
- The scale is heavy. Use caution, as lifting may cause it to fall over.

Consider the following installation conditions to get the greatest performance from your scale.

- The ideal conditions for installation are stable temperature and humidity, solid and level surface, location with no draft or vibration, indoors out of direct sunlight and a stable power supply.
- Do not install on a soft floor or where there is vibration.
- Do not install where there is wind or large fluctuations in temperature.
- Avoid locations in direct sunlight.
- Do not install in locations with strong magnetic fields or strong radio signals.
- Do not install where static electricity is likely to occur. When humidity is 45% R.H. or less, plastic and insulating materials are apt to be charged with static electricity due to friction, etc.
- Unstable AC power supply may cause malfunctions.
- Remove the protective film from the weighing pan before use.
- Sensitivity adjustment should be performed for more accurate weighing when the scale is first installed or when it is moved far. Refer to the [Instruction Manual: FG-CWP Series] (<https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html>) on our website <https://www.aandd.jp>.

2.2. Precautions for Storing

- Do not disassemble the scale.
- If the scale becomes dirty, wipe it with a soft cloth slightly moistened with a mild detergent or a rubbing alcohol (70% or less). Do not use organic solvents.
- Do not scrub with a brush or the like.
- Do not spray with strong jets of water.

2.3. Precautions for Weighing

- Do not place loads on the weighing pan that exceed the weighing capacity.
- Periodic sensitivity adjustment is recommended in order to maintain accurate weighing.
- Do not apply shock to or drop anything on the weighing pan.
- Do not use a sharp instrument such as a pencil or pen to press keys or switches.
- Press the **ZERO** key before each weighing to reduce the chance of error.
- Do not weigh with the scale immersed in water.
- Periodically confirm that weighing values are correct.
- The panel on the rear of the display unit should be kept closed when using the scale in order to make it dustproof and waterproof.


3. Items Included


The following items are packed.

- Main Unit


FG-30KCMWP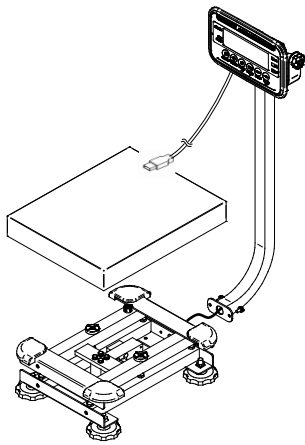

FG-60KCLWP
FG-150KCLWP

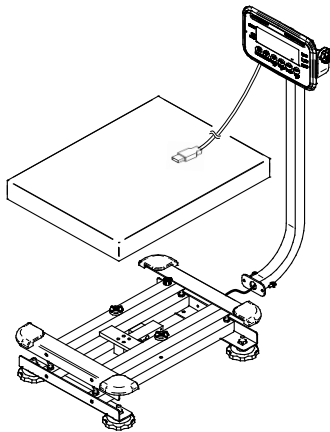- Quick Start Guide

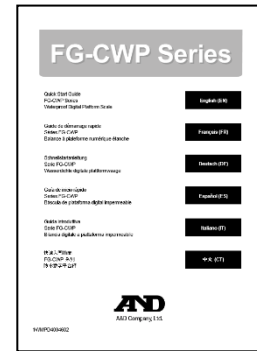
- AC Adapter

(Image diagram)


4. Part Names

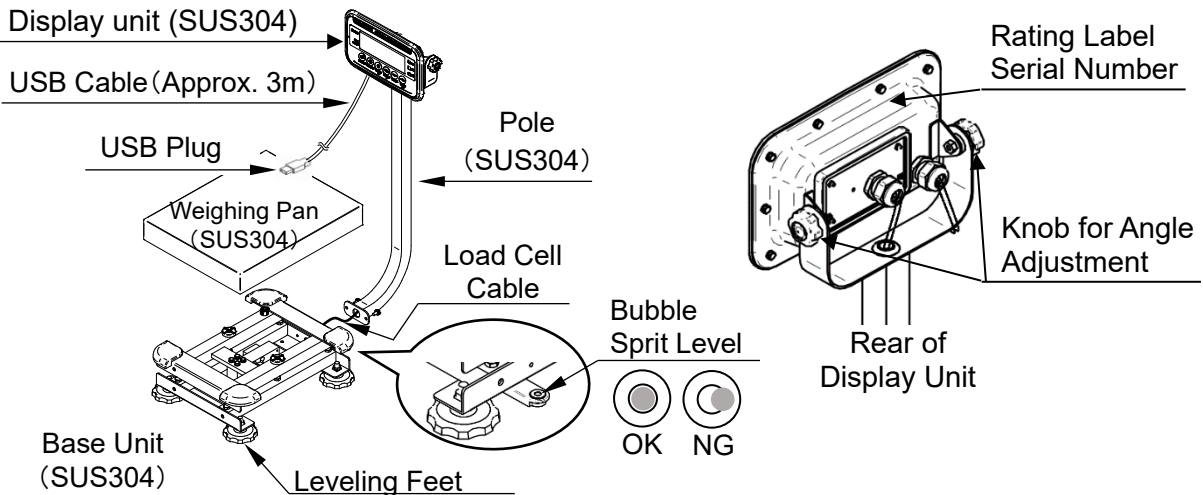4.1. Display and Symbols

Display Unit

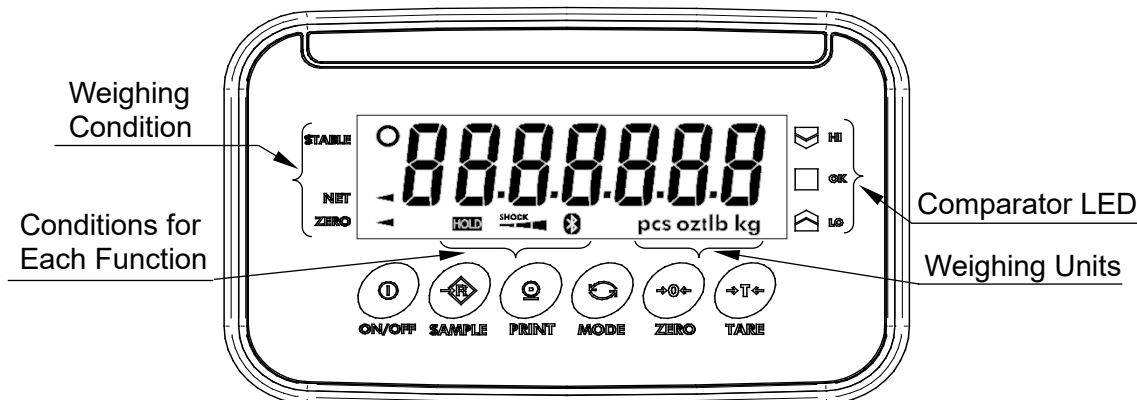
Display/symbol	Description
STABLE ○	This is lit when the weighing value is stable, indicating that the scale is in the proper condition for reading weighing values.
NET ◀	This is lit when weight of the container (tare) is subtracted.
ZERO ◀	This is lit when the scale is at the zero point (reference point for weighing).
Weighing Units	“kg”, “g”, “pcs”, “oz” and “lb” is lit. Note: “g” is only used with the FG-30KCMWP.
HOLD	This is lit when the display is held.
SHOCK	This is a function to detect impact to the mass sensor section and to display the impact level.
	Lights up when successfully connected to the optional FG-27CWP (Bluetooth communication interface).
HI OK LO	While the comparator function is being used, the weighing value is compared using the preset threshold values and the indicator displays the result.

- For details on the method of operation and various functions, refer to the [Instruction Manual: FG-CWP Series] (<https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html>) on our website <https://www.aandd.jp>.



4.2. Operation Keys

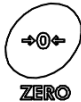Operation key	Description
	ON/OFF Key Shows or hides the display. Note: When the power plug of the scale is connected, power is supplied to the electric circuits inside the scale.
	SAMPLE Key Selecting "pcs" as the unit activates individual weight setting mode. Pressing and holding (for about 2 seconds) activates the comparator threshold setting mode
	PRINT Key Outputs the weighed value as data. Increases the value of the digit flashing when setting.
	MODE Key Switches the weighing unit. Shifts to the right the digit flashing when making settings.
	ZERO Key Based on the zero point weighed at the time of turning on the display, by pressing the ZERO key when the weight value is stable within $\pm 2\%$ of the weighing capacity, the zero point indicator turns on as soon as the display becomes zero.
	TARE Key The weight of the item on the weighing pan is subtracted as the weight of a container (tare weight).


5. Setup

1. Take the base unit and the pole out of the packing box while being careful not to pull the load cell cable.
2. Remove the weighing pan from the base.
3. Assembling the pole
 Temporarily remove the screws attached to the bottom of the weighing base, and then fasten the pole and base with the removed screws, taking care that the load cell cable is not pinched. Bundle and insert the excess load cell cable into the pole from the display side.
4. Put the weighing pan on the base unit.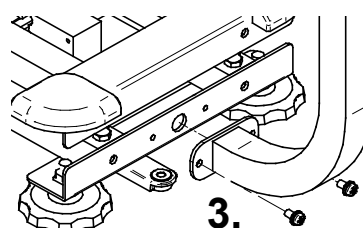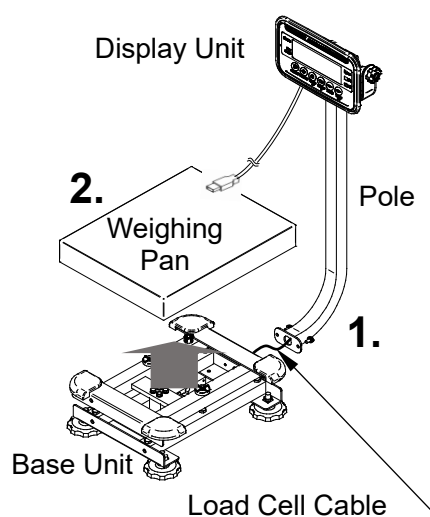

EN

- ⚠ When using in a place where static electricity is likely to be generated or in a wet place, join the grounding wire to the protective grounding terminal (screw).

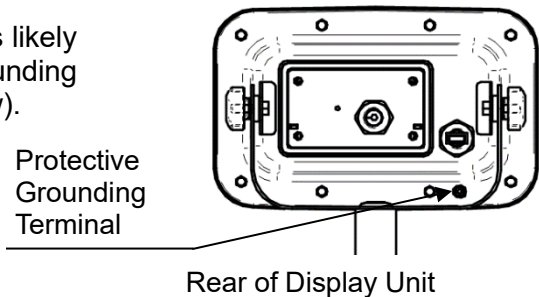6. Maintenance

- Refer to "2.Cautions" regarding use.
- Refer to "6.2.Error display" and corresponding mode for displayed error code.
- Refer to the [FG-CWP Series Instruction Manual] (<https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html>) our website <https://www.aandd.jp> for sensitivity adjustment.
- Periodically check the accuracy of weighing. Adjust the scale, if it is moved to another location or the environment has changed.

6.1. Repair

Do not disassemble / assemble the scale without an authorized service engineer. Doing so may cause an electric shock or damage to the scale, etc. In this case, repair is not covered under warranty. Contact your local A&D dealer if your scale needs service or repair.

6.2. Error display

In this situation	Confirm these items
Nothing displayed. Scale does not turn on.	<ul style="list-style-type: none"> • Is the main power the correct voltage? • Is the main power cord properly connected?
The scale does not display zero at first.	<ul style="list-style-type: none"> • Check around the weighing pan. • Is there anything on the weighing pan? • Refer to the [FG-CWP Series Instruction Manual] (https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html) on our website https://www.aandd.jp and perform zero point sensitivity adjustment.
E is displayed.	<ul style="list-style-type: none"> • Weighing error that meaning "Over loaded."
Lo ut is displayed.	<ul style="list-style-type: none"> • Indicates that the sample weight is too light to set the unit weight in the counting mode.
-CAL E is displayed.	<ul style="list-style-type: none"> • Sensitivity adjustment error that meaning "Too light."
Lb is displayed.	<ul style="list-style-type: none"> • Displayed when the main power drops.
----- is displayed and does not proceed.	<ul style="list-style-type: none"> • The weighed value is unstable due to drift, vibration or other factors. A breeze or vibration may be affecting measurement. Check around the weighing pan. • Check the connection of the load cell cable. • No zero display when the display is turned on. Remove anything that is on the pan. • Refer to the [FG-CWP Instruction Manual] (https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html) on our website (https://www.aandd.jp) and perform zero point sensitivity adjustment.

Fixed display.	<ul style="list-style-type: none"> • Did you use the "hold function" ? Refer to the [FG-CWP Series Instruction Manual] (https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html) on our website (https://www.aandd.jp) and change the function Hold. • Turn off scale and turn it on again.
----------------	---

6.3. Error code table

When any of the following errors are displayed, try turning the display off and on again.

Display	Content
Error0 is displayed	The temperature sensor has failed.
Error3 is displayed	The memory (circuit) has failed.
Error4 is displayed	The internal circuitry has failed.
Error5 is displayed	The mass sensor has failed.

7. Specifications

Models		FG-30KCMWP	FG-60KCLWP	FG-150KCLWP
kg	Weighing Capacity	30	60	150
	Readability	0.01	0.02	0.05
		0.005 *	0.01 *	0.02 *
		0.002	0.005	0.01
	0.001	0.002	0.005	
g	Weighing Capacity	30000	/	/
	Readability	10		
		5 *		
		2		
	1			
lb	Weighing Capacity	66	130	330
	Readability	0.02	0.05	0.1
		0.01 *	0.02 *	0.05 *
		0.005	0.01	0.02
	0.002	0.005	0.01	
oz	Weighing Capacity	1050	2,100	5,200
	Readability	0.5	1	2
		0.2 *	0.5 *	1 *
		0.1	0.2	0.5
	0.05	0.1	0.2	
lb-oz	Weighing Capacity	66	130	330
	Readability	0.1	1	1
Number of Samples		5 pcs (can be changed to 10 pcs , 20 pcs, 50 pcs or 100 pcs)		
Maximum Count Number		300,000 pcs		
Minimum Unit Mass [kg]		0.0001	0.0002	0.0005
Repeatability (Standard Deviation) [kg]		0.005	0.01	0.02
Linearity [kg]		±0.01	±0.02	±0.05
Stabilization time (good environment)		20kg:Approx. 1.0 sec (Factory setting) Approx. 0.8 sec (Land 0 setting)		

Models	FG-30KCMWP	FG-60KCLWP	FG-150KCLWP
Temperature Drift	±20 ppm/°C typ. (5 °C to 35 °C)		
Display	7 segment LCD with backlight(Character height 30 mm)		
Display update	Approx. 10 times/second		
Dust and water protection	Complying with IP67		
Operating Conditions	-10 °C to 40 °C, 85% R.H. or less (non-condensing)		
Power Supply	AC adapter or mobile battery or USB port USB Cable Length: Approx. 3 m (USB cable is also used for communication.)		
Weighing Pan Size [mm]	300 x 380	370 x 500	
Dimension [mm] Width x Depth x Height	300 x 515 x 784	370 x 635 x 784	
Weight (approximately)	11.1 kg 24.5 lb	12.8 kg 28.3 lb	
Sensitivity adjustment weight (factory setting)	30 kg 60 lb	60 kg 120 lb	150 kg 300 lb

* **Factory setting**

FG-CWP Series Class III Model Specifications

Models		FG-30KCMWP	FG-60KCLWP	FG-150KCLWP
Accuracy class		III		
kg	Maximum Capacity	30	60	150
	Readability	0.01	0.02	0.05
lb	Maximum Capacity	66	130	330
	Readability	0.02	0.05	0.1
oz	Maximum Capacity	1,050	2,100	5,200
	Readability	0.5	1	2
Minimum Unit Mass [kg]		0.2	0.4	1
Maximum tare [kg]		30	60	150

- Class III approved model specifications. For details, refer to the [Instruction Manual: FG-CWP Series] (<https://www.aandd.jp/products/weighing/scale/platform/fgcwp.html>) on our website <https://www.aandd.jp>. The appearance is the same as that of the FG-CWP series.

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A&D AUSTRALASIA PTY LTD
A&D KOREA Limited
A&D Instruments India Private Limited