HL-i series Digital Compact Scale Operating Instructions

1WMPD4000944C

Cautions

- Avoid installing the scale in direct sunlight.
- Avoid over / shock load that could cause damage to the scale.
- Use the scale under conditions of -10°C to 40°C, less than 85% R.H.
- The weighing pan cannot be removed. Do not pull up on it.
- Do not mix the battery types. Replace all of the batteries at same time.
- If the scale is not to be used for a long period of time, remove the all batteries.
- Use only the specified AC adapter for the scale (AC adapter is optional).

Compliance with FCC Rules

Please note that this equipment generates, uses and can radiate radio frequency energy. This equipment has been tested and has been found to comply with the limits of Class A digital devices pursuant to Part 15 of FCC rules. These rules are designed to provide reasonable protection against interference when equipment is operated in a commercial environment. If this unit is operated in a residential area, it may cause some interference and under these circumstances the user would be required to take, at his own expense, whatever measures are necessary to eliminate the interference. (FCC = Federal Communications Commission in the U.S.A.)

Specifications

| | |
|--------------|---|
| HL-200i | 200 x 0.1 g / 7.05 x 0.01 oz / 1000.0 x 0.5 ct / 5.335 x 0.005 tl |
| HL-2000i | 2000 x 1 g / 70.5 x 0.1 oz / 4 lb 6.6 oz x 0.1 oz / 53.35 x 0.05 tl |
| Power | 4 x "AA" size batteries (monitor batteries included) or AC adapter |
| Battery life | Approximately 1000 hours with alkaline cells at 20 °C / 68 °F |
| Pan size | φ 130 mm / φ 5.1 in. |
| Dimensions | 130 (W) x 192 (D) x 51 (H) mm / 5.1 (W) x 7.6 (D) x 2.0 (H) in |


Keys and Symbols

| | |
|-----------------|---|
| | Turns the scale power on/off. |
| | Sets the display to zero. |
| | Changes weighing units. |
| Symbol STABLE ● | Indicates when the reading is stable. |
| Symbol ZERO ◀ | Indicates when the scale zero is correct. |

Installing batteries / Connecting the AC adapter

Installing / Replacing Batteries

Remove the battery cover and insert 4 batteries (AA size) into the battery housing, especially take care of the polarities. Attach the battery cover again.

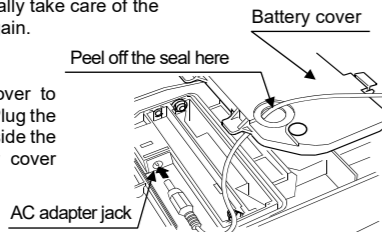AC adapter (optional)

Peel off the seal on the battery cover to thread the AC adapter wire through. Plug the AC adapter to the AC adapter jack inside the battery housing. Attach the battery cover again.

Weighing procedures

1. Turn on / off the scale

Press the key to turn on the scale. The scale displays all segments for a few seconds and then displays zero. Press the key again to turn the scale off.

2. Select the weighing unit

Press the key to select a weighing unit.
HL-200i: "g" → "oz" → "ct" → "tl" → "g" "tl" = Taiwan tael
HL-2000i: "g" → "oz" → "lb-oz" → "tl" → "g"

3. Weighing

Verify the reading is zero. If not, press the key to zero the display. If you use a container for weighing, place an empty container on the weighing pan and press the key to zero the display. Place the objects to be weighed on the pan or in the container. When the reading becomes stable, the STABLE indicator ● is displayed.

Note

Automatic power-off function

If the scale is left ON and the STABLE indicator ● is displayed, the automatic power-off function turns the scale off after approximately 5 minutes. To disable or enable this function, press the key while pressing the key. When the scale is turned on, (XXX = 1.00 etc.) is displayed and this function is disabled. If is not displayed, then the function is enabled.

Error message

| | |
|--|---|
| | Overload. Remove the objects from the pan. |
| | Low battery. Replace the batteries with new ones. |

Calibration

For information about calibration, please visit the A&D website (<https://www.aandd.jp>).

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