

# Connecting Viscometers to the AD-8127 Compact Printer

2017.07.20

## 1. Introduction

This document describes usage and setting examples for connecting viscometers and the AD-8127 compact printer. Refer to the respective manual for how to change settings of the viscometers and the printer.

## 2. Applicable models

Viscometers

SV-10, SV-100, SV-1A, SV-10A, SV-100A

Rheometers

RV-10000A, RV-10000

## 3. Usage and setting examples

### Usage example

Printing the measurement results.

Viscometer internal settings			AD-8127 Compact printer Printing mode
Data output mode Prt	Data output format tYPE	Measurement elapsed time output S-At	
0, 1	1	0, 1	DUMP
Date/time output S-td	Other output S-Ed	Pause at data output PUSE	
0, 1	0, 1, 2	0	