


How to connect to the balance


AD-8922

Remote Controller

- Applicable balances are listed in the table below.
 ※ For other models (setting procedure/key operation), refer to the relevant manual.
- Connect the balance to the AD-8922, using the communications cable provided with the AD-8922, and turn them both on.
- Set the balance data output mode to "stream mode" (Continuous output).
 ● For GX,GX-K,GF,GF-K,GP,GH,GR,HR-i,EK-H, set the function setting to "dout Prt 3".

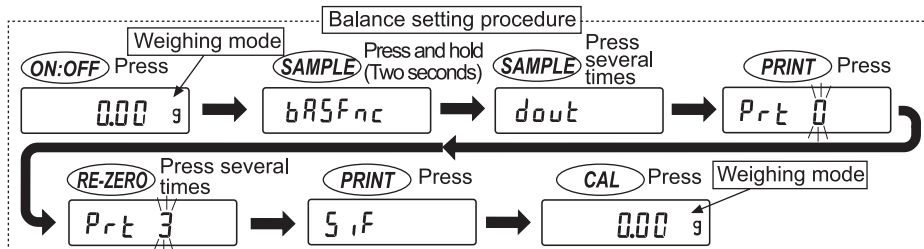

- Referring to the table below, change the AD-8922 function setting of "cSEt".

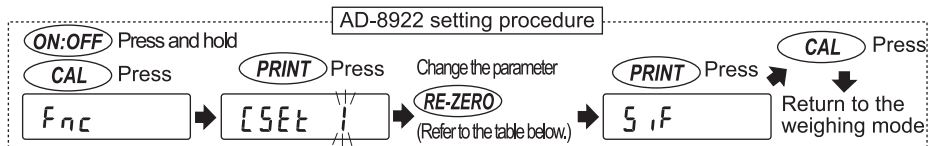

- AD-8922 key operation (Depends on the balance)

Weighing instrument	Command set	AD-8922 key					
		ON:OFF	CAL	SAMPLE	PRINT	MODE	RE-ZERO
GX,GX-K,GP,GH	cSEt 1	Turns the weighing instrument display on or off.	Calibrates using the internal mass.	Switches the minimum display.	Outputs the received data to an external device.	Switches the unit displayed.	Sets the display to zero.
GR	cSEt 2		—				
GF,GF-K,EK-H,HR,HR-i	cSEt 3	—	—	—	—	—	—
EK-i,EW-i	cSEt 4	—	—	—	—	—	—