

Additional Explanation for the AD-1690

This document is additional information for the instruction manual of the leak tester AD-1690. Make use of this additional explanation in addition to the AD-1690 instruction manual.


Additional contents

When your hand tip can not be attached to the attachment due to its special shape, use the adapter tube.

The following explains how to use this adapter tube.

About accessories

With the “Adapter and Attachment” in the “3.2. Accessories” on page 4 in the AD-1690 instruction manual, the adapter tube is added and the constitution is as follows.

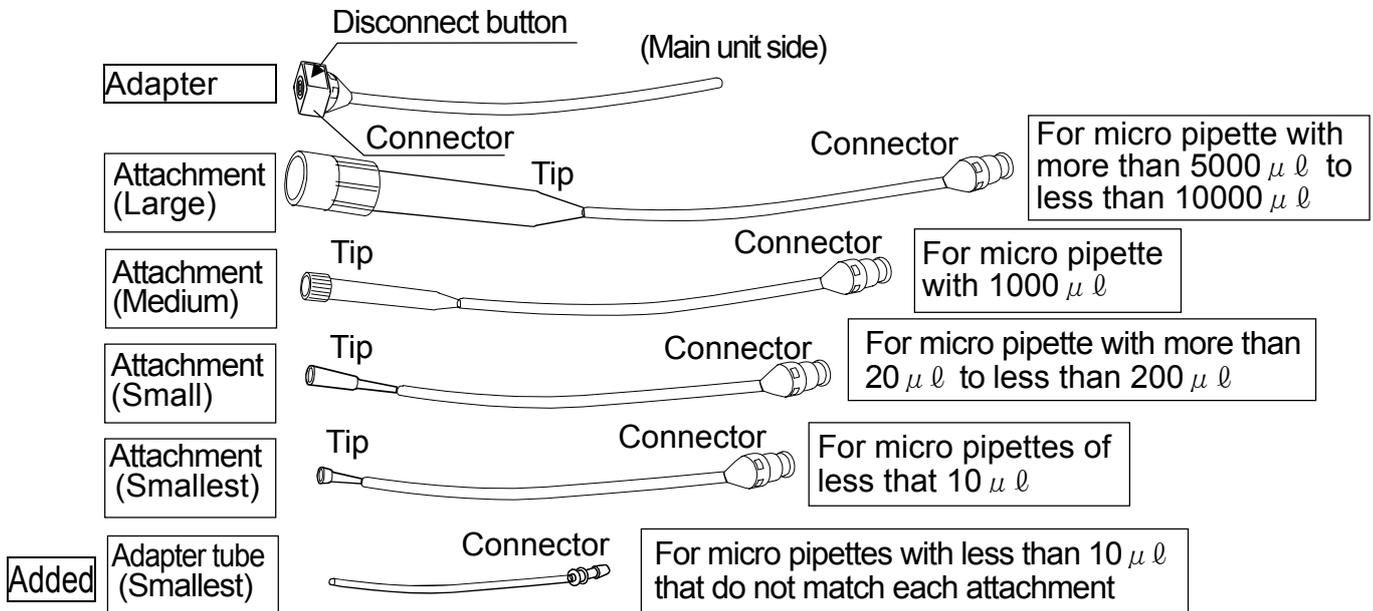

Using the adapter tube

By following the “(2) Connecting the attachment (pipette side)” in the “5. MEASUREMENT PREPARATION” on page 7, remove the tip from the provided attachment and exchange with your special tip.

