

WinCT (Windows Communication Tools) Ver. 4.01 for Windows Vista
English edition
CD-ROM for windows
Copyright (c) 1998 A&D Company limited.

Setup text

WinCT (Windows Communication Tools)

WinCT is a program for transmitting data from an A&D balance to a computer.
Below are instructions for installation and use of WinCT.

1. System Requirements

Computer OS Microsoft Windows Vista / XP
 RS-232C 1port
 - If the computer has no COM port, a Serial/USB converter is
 available as an accessory to add a COM port on the computer.
 A cable will also be attached for the connection with a balance.
 The balance's RS-232C has 25 pins -> Model No. AX-USB-25P-EX
 The balance's RS-232C has 9 pins -> Model No. AX-USB-9P-EX

Balance An A&D balance with an RS-232C interface
Cable An RS-232C cable

CAUTION

- Use an English version of Windows for proper operation of WinCT.
- If WinCT is operated on Windows except for an English version, the computer may show incorrect characters of data which is already stored before installing WinCT.
- Back up your important data in the computer before installing WinCT.

- For Windows XP, Microsoft .Net Framework 2.0 or higher has to be installed. (Microsoft .Net Framework is a freeware available on Microsoft's website.) If Microsoft .Net Framework is already installed, when you open "Add or Remove Programs" in the "Control Panel," it will appear in the "Currently Installed Programs" list.

2. Installation Instructions

- Step 1 If a previous version of WinCT is installed on the computer, uninstall the previous version of WinCT.
- Step 2 Close all open applications in Windows.
- Step 3 Install WinCT.
 Run WinCT_Setup.msi.
- Step 4 Install WinCT following the instructions on the screen.

3. Starting WinCT

Select either RsCom or RsKey from the START menu. The default menu address is [Start] - [program] - [A&D WinCT]. You will find the instruction manual in the Readme file under [A&D WinCT].

RsCom This window is able to transmit the weighing date using a text file format to a computer. This window also transmits commands for operating an A&D balance.

RsKey Weighing data is imported directly into an application, such as MS Excel. However, this window cannot be used to send commands to a balance.

CAUTION

- WinCT is a program for balances that use the A&D format. A&D does not guarantee performance or compatibility when other manufacture's balances are used with WinCT.
- If you have problems or questions about compatibility with third party software, contact the manufactures for information.

- This version does not include the real-time graph plotting software, "RsWeight."

4. License Agreement and Limited Warranty

- WinCT is subject to change without notice and does not represent a commitment by A&D to support earlier versions or to make changes to support non standard computers.
- WinCT is intended for the transmission of the weighing data from an A&D balance to a PC, and command transmission for controlling the balance.
- Under copyright laws, the WinCT program may not be copied, in whole or part, without the consent of A&D, except as specified in the license agreement.
- WinCT should only be installed on the hard disk or peripheral storage devices of a computer connected to an A&D balance.
- A&D bears no liability for direct, indirect, special, incidental, or consequential damages resulting from any defect in the WinCT program or the readme file, even when advised of the possibility of such damage.
- A&D is not responsible or liable for the loss of any programs and/or data stored in the computer on which WinCT was installed, including the costs of recovering said lost programs or data.

Microsoft, Windows, Excel and Word are trademarks or registered trademarks of Microsoft Corporation.

5. Uninstall Instructions

Use the "Add/Remove Programs" icon in the control panel to remove this program from your computer.

End of Setup Text
