

Verified Operation of Our Software & Devices for Weighing Products with Windows 7 (32-bit Version)

Updated: January 2010

Description		Results
WinCT Ver.3.00	RsCom	Verified
	RsKey	Failedⁱ
	RsWeight	Verified
WinCT Ver.4.01	RsCom	Verified
	RsKey	Verified ⁱⁱ
	RsWeight	Not available ⁱⁱⁱ
WinCT-Moisture Ver.2.41	RsFig	Verified
	RsTemp	Verified
WinCT-Viscosity Ver.1.11	RsVisco	Verified
WinCT-Plus Ver.2.10	RsMulti	Verified
	DeviceInstaller	Verified ^{iv}
	Com Redirector	Verified
WinCT-AD4212C Ver.1.02		Verified
WinCT-Pipette Ver.1.1.4		Verified
AD-1688		Verified
AD-8527		Verified
OP-02 (Quick USB)		Verified
AX-USB-xx (Serial/USB converter)		Verified ^v

ⁱ Data cannot be transferred.

ⁱⁱ The data transfer rate was slower compared with Windows Vista/XP.

Windows Vista/XP: Approx. 6.3 times/sec.

Windows 7: Approx. 1.4 times/sec.

ⁱⁱⁱ Please use the version 3.00 should this function be required.

^{iv} As with Windows Vista, Excel cannot be started from RsMulti.

^v Please use the driver software for Windows Vista.