


1. Features

The FX-10 protects the balance from the effects of a breeze and will provide more accurate weighing results.

All components are antistatic, making it less likely that weighing errors will occur due to static electricity.

Models : FZ / FX, FZ-WP / FX-WP


2. Unpacking

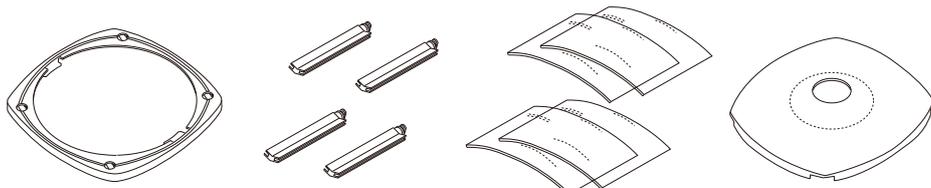

Base 1 pcs Pillars 4 pcs Plates 4 pcs Cover 1 pcs



3. Assembling

1. Insert the pillars into the holes in the base.
2. Rotate the pillars 90 degrees, and confirm that the position of the channels on the pillars and the base match up.
3. Insert the plates in the channels between the pillars and aligned with the base.
4. Confirm that the projections on both sides of the bases are aligned with the channels on both sides of the balance. Install the breeze break on the balance.
5. Place the cover on the breeze break.

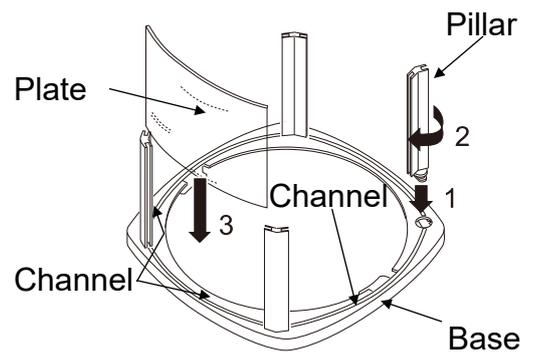

Note (FZ-WP / FX-WP series)

When cleaning the balance, remove the small breeze break.

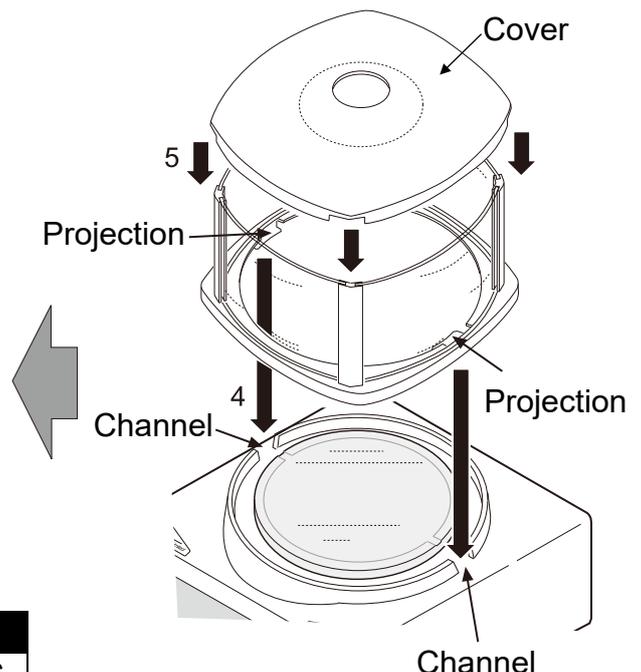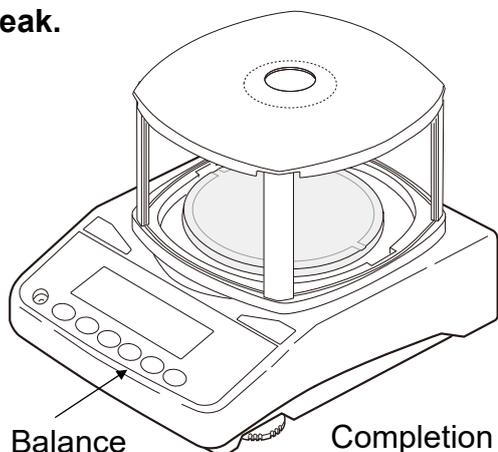

Materials used (Material of the breeze break)
 Antistatic treated Acrylic (PMMA), Antistatic ABS

* Each part has the type of material used printed on it.