

GP-12


Animal Weighing Bowl

Instruction Manual

Models

GX-L/GF-L series

GP series

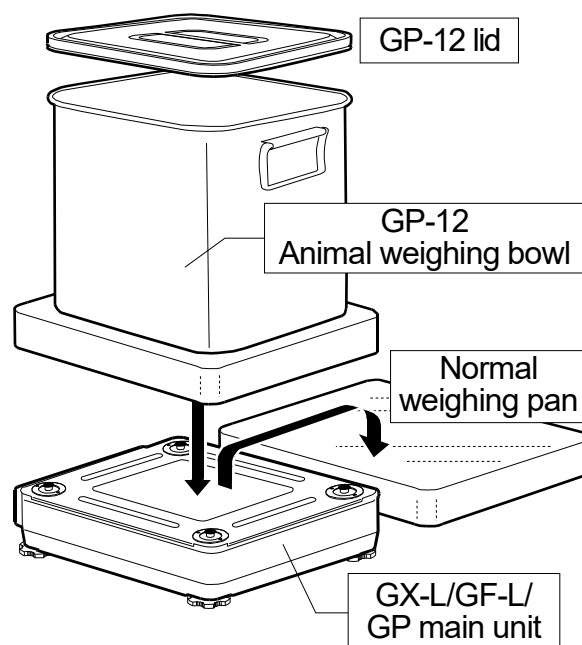

Weighing

Setup

- 1 Replace the normal weighing pan with the animal weighing bowl (GP-12).

Measurement

- 2 • GX-L/GF-L series:
Press the **ZERO** key or the **TARE** key to cancel the weight (tare).
• GP series:
Press the **RE-ZERO** key to cancel the weight (tare).
- 3 Place a sample animal in the bowl to weigh it.

Caution

- When using the animal weighing mode (Hold function), the average weight of the animal can be displayed. For details, please refer to the instruction manual.
- When using this option, the weighing range that can be used is an approximately 4 kg less than the weighing capacity.
- If necessary, use the GP-12 lid.
- When performing sensitivity adjustment on the balance, use the normal weighing pan.
- When performing sensitivity adjustment on the balance with the animal weighing bowl and replacing with the normal weighing pan, **-E** is displayed. If the normal weighing pan is used, re-adjust the balance sensitivity with it.