


GXK-012

Animal Weighing Bowl

Instruction Manual

Models

GX-K series
GF-K series

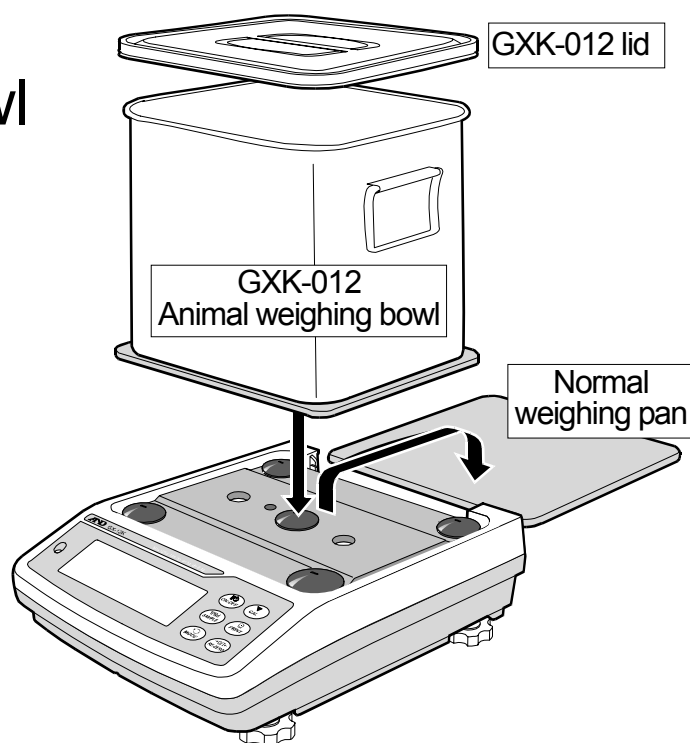

Weighing

Setup

- 1 Replace the normal weighing pan with the animal weighing bowl (GXK-012).

Measurement

- 2 Press the **RE-ZERO** key to cancel the weight (tare).
- 3 Place a sample animal in the bowl to weigh it.

Caution

- When using the animal weighing mode (Hold function), the average weight of the animal can be displayed.
- When using this option, the weighing range that can be used is an approximately 1.5 kg less than the weighing capacity.
- If necessary, use the GXK-012 lid.
- When calibrating the balance, use the normal weighing pan.
- When calibrating the balance with the animal weighing bowl and replacing with the normal weighing pan, **-E** is displayed. If the normal weighing pan is used, re-calibrate the balance with it.