


GXK-012

Animal Weighing Bowl

Instruction Manual

Models

GX-K series
GF-K series
GX-M series
GF-M series

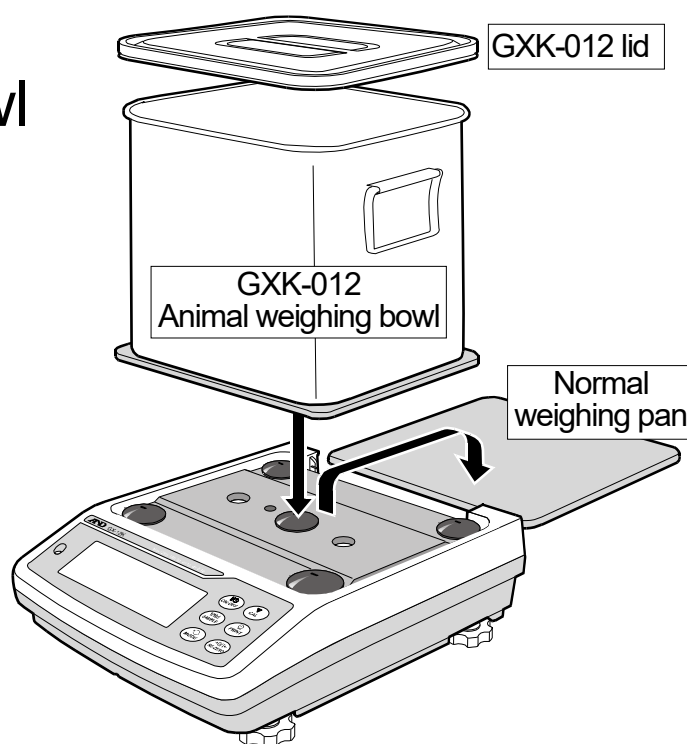

Weighing

Setup

1. Replace the normal weighing pan with the animal weighing bowl (GXK-012).

Measurement

2. Press the **RE-ZERO** key to cancel the weight (tare).
3. Place a sample animal in the bowl to weigh it.

Caution

- When using the animal weighing mode (Hold function), the average weight of the animal can be displayed.
- When using this option, the weighing range that can be used is an approximately 1.5 kg less than the weighing capacity.
- If necessary, use the GXK-012 lid.
- When performing sensitivity adjustment with the balance, use the normal weighing pan.
- If performing sensitivity adjustment with the balance that has the animal weighing bowl on it and then replacing it with the normal weighing pan, **-E** is displayed. When using the normal weighing pan again, re-perform sensitivity adjustment with the balance that has the normal weighing pan on it.

AND

A&D Company, Ltd.