
Compact Scale : HL-100

INSTRUCTION MANUAL

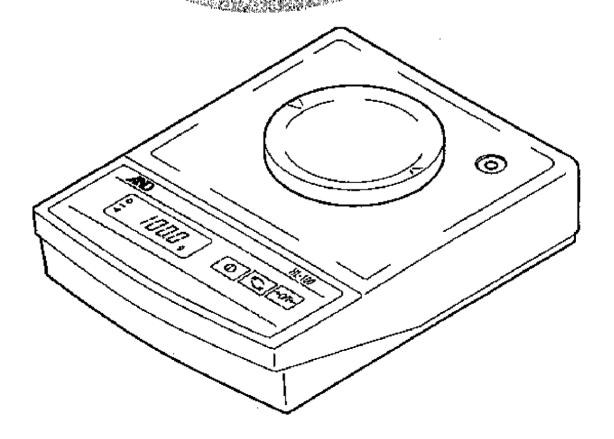



SAFETY PRECAUTIONS

All safety messages are identified by the following the words " WARNING " and "CAUTION". These words mean the following:

⚠ WARNING	Important information to alert you to a situation that might cause serious injury and damage to your property if instructions are not followed.
⚠ CAUTION	Important information that tells how to prevent damage to equipment.

When using the HL, the following safety precautions should always be followed.

Λ WARNING

Use only the AC adapter which comes with the scale. Other adapter may cause damage. (AC adaptor is optional)

Internal service or adjustment to this product should be performed by a qualified person.

\triangle CAUTION

Avoid installing the scale in direct sunlight, that may cause discoloration or malfunctions.

Do not mix the type of batteries. Replace all other batteries at the same time.


If the scale is not to be used for a long period of time, remove all batteries from the battery compartment to avoid leakage which may cause damage to the instrument.

Avoid overloading that could cause damage to the scale.

Do not drop water in the scale which is not water-resistant. It causes damage, if inside of the scale is wet.

The weighing could be influenced by matter charged static electricity. Discharge the static electricity. As example, the method is to use the electrification prevention spray, to spray it to both side of weighing platform.


PARTS DESCRIPTION

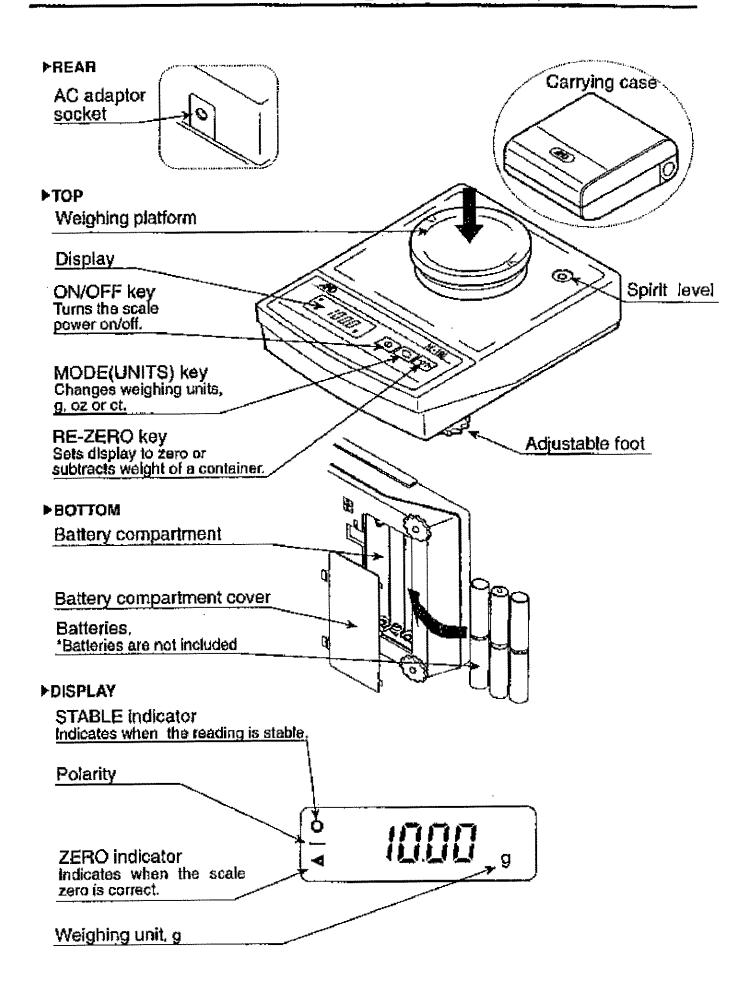

PREPARATION


1. Installing batteries / Connecting the AC adaptor

Batteries

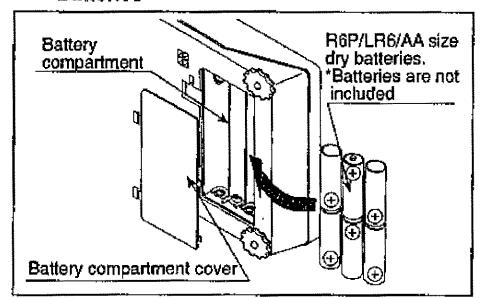

Remove the battery compartment cover and insertsix batteries (R6P/LR6/AA size) into the battery compartment, taking extreme care that the polarities plus and minus are observed.

AC adaptor (optional)

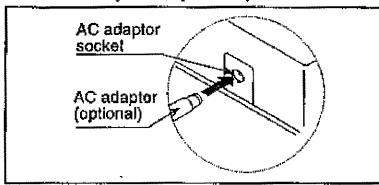

Plug the AC adaptor to the AC adaptor socket on the side. The AC input requirement could be 100, 120, 220,230, or 240Volts (50/60Hz) depending on the area where used, so please verify that the adaptor is correct.

2. Setting up the HL

A CAUTION

Avoid installing the scale in direct sunlight, that may cause discoloration or malfunctions. Place your HL on a firm weighing table so that the scale is level. (The scale will not perform accurately when it is not level.)

Place the HL on the firm surface and adjust the feets so that the spirit level shows the scale is level for an accurate weighing.

WEIGHING

Before weighing

Be aware of the Automatic power-off function.

If this unit is left ON and the stable indicator is displayed, an automatic power-off function turns power off after approximately 5 minutes. To disable this function, press [ON/OFF] while pressing [RE-ZERO]. "P-XX" will be displayed and the scale returns to the weighing mode.

Error messages

F Overload

---- The scale zero is out of range.

Lb Low battery

Calibration may be required before weighing.

Read "CALIBRATION" first and if necessary, calibrate your scale for accurate weighing.

Weighing procedures

1. Press [ON/OFF] to turn on the scale.

When power is turned ON, all display segments appear for a few seconds and "0" will appear on the display.

2. Select the weighing unit with [MODE] ([UNITS]).

Press [MODE] ([UNITS]) to select a weighing unit either "g", "oz" or "ct".

Once the unit has been selected, the selected unit will be displayed to the right of the weight value.

3. Start weighing

If you do not use a container for weighing,

Verify the reading is "0". If not, press [RE-ZERO] to display "0".

Place objects on the weighing platform to weigh.

When the reading becomes stable, the stable indicator is displayed.

If you use a container for weighing,

Place an empty container on the platform.

Wait for the stability indicator to be displayed and press [RE-ZERO].

Place the objects to be weighed in the container.

When the reading becomes stable, the stable indicator is displayed.

Select the weighing unit

HL-100 scale has "g", "oz" or "ct" selection. Scale is shipped from the factory to power up in "g". To select "oz" or "ct" mode, turn scale off. Press [ON/OFF] While pressing [MODE]. Scale will power up in "oz"

or "ct" mode. Repeat this instruction, then power-up unit changes cyclically "g" \Rightarrow "oz" \Rightarrow "ct" \Rightarrow "g".

Note) "oz" is available in non-metric countries only.

CALIBRATION

When to calibrate

Calibration may be required when it is initially installed, if the scale is moved a substantial distance, or in accordance with local regulations. This is necessary because the weight of a mass in one location is not necessarily the same in another location. Also, with time and use, mechanical deviations can occur.

Calibration using a calibration weight

Calibrate a scale with calibration weights at capacity or 1/2 capacity. 100g \pm 0.002g ("CAL F") 50g \pm 0.002g ("CAL5")

1. Enter calibration mode.

Press [ON/OFF] to turn the power off. While pressing both [RE-ZERO] and [MODE] ([UNITS]), press [ON/OFF]. "CAL" will be displayed.

2. Zero calibration

Press [RE-ZERO] to calibrate zero. "CAL 0" will be displayed. Wait for the stable indicator to be displayed and press [RE-ZERO] "CAL F" will be displayed in a few seconds.

3. Span calibration

To calibrate by using the 50g weight, in the "CALF" display, press the [MODE] ([UNITS]) to display "CAL 5".

Place the correct weight at the center of the platform for an accurate calibration.

Wait for the stable indicator to be displayed, then press [RE-ZERO]. The display will show "END" and automatically the scale return to the weighing mode.

Calibration by gravity compensation

If you have calibration weights at the capacity of the scale, you can calibrate the scale with the weights instead of compensating for acceleration due to gravity.

The HL is calibrated at 9.798 m/sec². If the acceleration of gravity at your location is not the same, calibrate the scale by compensating the

acceleration of gravity. (See page 7 "The Value of Gravity at Various Locations" and page 8 "World Map" and find your location' value.)

1. Enter the calibration mode.

Press [ON/OFF] to turn the power off. While pressing both [RE-ZERO] and [MODE] ([UNITS]), press [ON/OFF]. "CAL" will be displayed.

2. Select a digit to be changed.

Press [MODE] ([UNITS]). "9.798" will be displayed. Press [RE-ZERO], the decimal point will move and the last digit will be selected. Select a digit to be changed with [MODE] ([UNITS]). The first digit after decimal point is the selected digit.

3. Set value.

Set the value of the digit selected by pressing [RE-ZERO]. (See "The Value of Gravity at Various Locations" and "World Map" and find your location' value.)

4. Enter the value in memory.

While pressing [MODE] ([UNITS]), press [RE-ZERO] to enter the value.

END will be displayed.

SPECIFICATIONS


MODEL	HL-100
Capacity	100 g / 3.5 oz* / 500 ct
Resolution	0.01 g / 0.0005 oz* / 0.05 ct
Non-linearity	±0.03 g
Repeatability	±0.02 g
Calibration weight	100 g or 50 g ±0.002 g
Span drift	±0.015%/°C TYP (5°C~35°C /41°F~86°F)
Display	8 mm/0.31 inch, 7segment liquid crystal display
Platform size	₱ 70 mm / ₱ 2.75 inches
Dimensions	140(W)×150(D)×55(H) mm, 5.5(W)×5.9(D)×2.17(H) inches
Weight	Approximately 460 g/1lb
Power	6 × R6P/LR6/ "AA" size batteries or AC adaptor
Battery life ·	Approximately 70 hours with manganese type
cells	/ 140 hours with alkaline cells at 20°C/68°F
Operating temp.	-10°C~40°C /14°F~104°F, Less than 85%RH (Non-condensing)
Accessories	This manual, Carrying case
Options	AC Adaptor

^{* &}quot;oz" is only for non-metric countries.


THE VALUE OF GRAVITY AT VARIOUS LOCATIONS

Amsterdam	9.813	m/s²
Athens	9.800	m/s²
Auckland NZ	9,799	m/s²
Bangkok	9.783	π/s²
Birmingham	9.813	m/s²
Brussels	9.811	m/s²
Buenos Aires	9.797	π/s²
Calculta	9.788	m/s²
CapeTown	9.796	π/s²
Chicago	9.803	
Copenhagen	9.815	m/s²
Cyprus	9.797	m/s²
Djakarta	9.781	m/s²
Frankfurt	9.810	m/s²
Glasgow	9.816	m/s²
H a vana	9,788	m/\$²
Helslnki	9,819	m/s²
Kuwait	9.793	m/s²
Lisbon	9.801	m/s²
London (Greenwich)	9.812	π/s²
Los Angeles	9.79 6	
Madrid	9.800	m/s²

Manila	9.784	rn/s²
Melbourne	9.800	m/s²
Mexico City	9.779	rn/s²
Milan	9.806	rrv/s²
New York	9.802	
Osio	9.819	
Ottawa	9.806	
Paris	9.809	m/s²
Rio de Janeiro	9.788	-
Rome	9.803	
San Francisco	9.800	
Singapore	9.781	
Stockholm	9.818	
Sydney	9.797	
Talchung	9.789	
Talwan	9.768	
Taipel	9.790	
Tokyo	9.798	m/s²
Vancouver, BC	9.809	
Washington DC	9.801	
Wellington NZ	9.803	
Zurich	9.807	

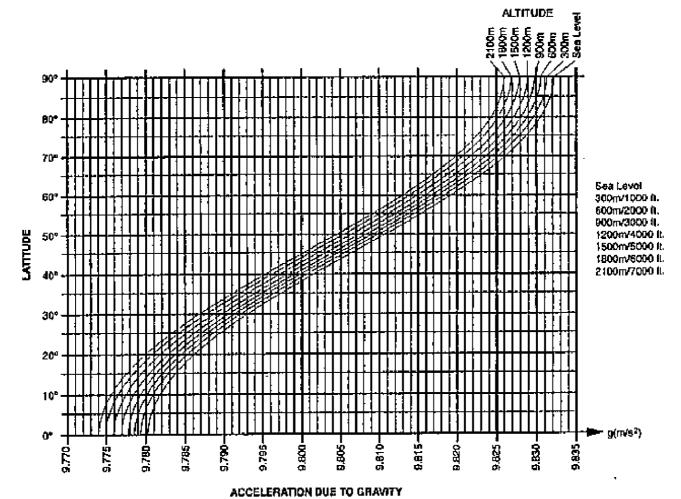

WORLD MAP

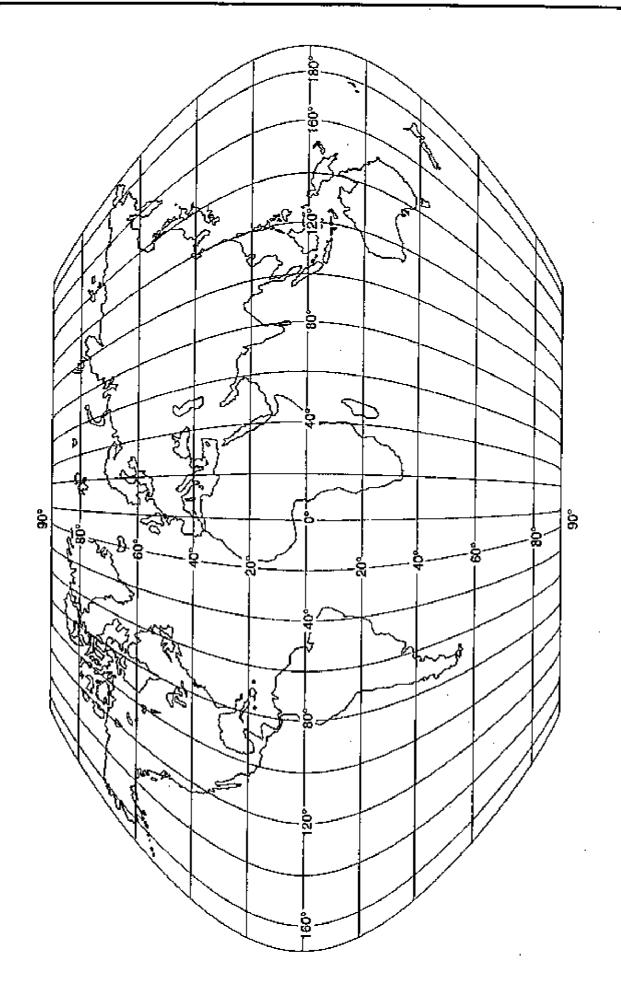



A&D Company, Limited

3-23-14 Higashi-ikebukuro, Toshima-ku, Tokyo 170-0013 JAPAN Telephone: [81] (3) 5391-6132 Fax: [81] (3) 5391-6148

A&D ENGINEERING, INC.

1555, McCandless Drive, Milpitas, CA. 95035 U.S.A. Telephone: [1] (408) 263-5333 Fax: [1] (408)263-0119

A&D INSTRUMENTS LTD.

Unit 24/26 Biacklands Way, Abingdon Business Park, Abingdon, Oxon OX14 1DY United Kingdom Telephone: [44] (1235) 550420 Fax: [44] (1235) 550485

<German Scales Office>

Berner Strabe 64, 60437 Frankfurt/Main 50 GERMANY Telephone: [49] (69) 507-1017 Fax:[49] (69) 507-2054

A&D MERCURY PTY. LTD.

32 Dew Street, Thebarton, South Australia 5031 AUSTRALIA Telephone: [61] (8) 8352-3033 Fax: [61] (8) 8352-7409

A&D KOREA Limited

8th Floor, Manhattan Bldg. 36-2 Yoido-dong, Youngdeungpo-ku, Seoul, KOREA Telephone: [82] (2) 780-4101 Fax: [82] (2) 782-4280