HT series Compact Scale Operating Instructions

1WMPD4001904D

Cautions

- Avoid installing the scale in direct sunlight.
- Avoid over / shock load that could cause damage to the scale.
- Use the scale under conditions of -10 °C to 40 °C, less than 85% R.H.
- Do not mix the type of batteries. Replace the batteries all together.
- If the scale is not to be used for a long period of time, remove all batteries.
- Use only the specified AC adapter for the scale (AC adapter is optional).

Compliance with FCC Rules

Please note that this equipment generates, uses and can radiate radio frequency energy. This equipment has been tested and has been found to comply with the limits of Class A digital devices pursuant to Part 15 of FCC rules. These rules are designed to provide reasonable protection against interference when equipment is operated in a commercial environment. If this unit is operated in a residential area, it may cause some interference and under these circumstances the user would be required to take, at his own expense, whatever measures are necessary to eliminate the interference. (FCC = Federal Communications Commission in the U.S.A.)

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Keys and Symbols

[ON/OFF] key	Turns the scale power on/off.
[RE-ZERO] key	Sets the display to zero.
[UNITS] key	Changes weighing units, g, oz or lb-oz.
Symbol	Indicates when the reading is stable.
Symbol	Indicates when the scale zero is correct.
Symbol NET	Indicates when the net weight is displayed.
Symbol	Battery remaining level, shown after power-on only.

Installing the batteries / Connecting the AC adapter

Installing / Replacing the batteries

Remove the battery cover.

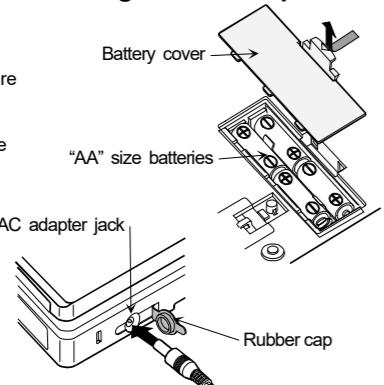
Insert 4 new batteries (AA size) into the battery compartment, taking extreme care of the polarities.

Replace the battery cover.

The batteries provided with the scale are for testing the performance of the scale and may have a limited life.

AC adapter (optional)

Verify that the AC adapter is correct. Open the rubber cap and plug the AC adapter into the AC adapter jack.

Weighing procedures

1. Turning the scale on / off

Press the [ON/OFF] key to turn the scale on. The scale displays all segments for a few seconds and then displays zero. Press the [ON/OFF] key again to turn the scale off.

2. Selecting the weighing unit

Press the [UNITS] key to select the weighing unit. The scale will power up with the weighing unit that was in use when turned off last time.

3. Weighing

Verify that the reading is zero. If not, press the [RE-ZERO] key to zero the display. If you use a container for weighing, place an empty container on the weighing pan and press the [RE-ZERO] key to zero the display. Place the objects to be weighed on the pan or in the container. When the reading becomes stable, the STABLE indicator is displayed.

Note

Automatic power-off function

If the scale is left ON and the STABLE indicator is displayed, the automatic power-off function turns power off after approximately 5 minutes. To disable or enable this function, press the [ON/OFF] key while pressing the [RE-ZERO] key.

Then

(disabled) or

Error message

Overload. Remove the objects from the pan.

The scale is out of zero range or unstable when powered on.

Low battery. Replace the batteries or use the AC adapter.

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