


Charging stand for four pipettes

AX-ST-CH-M4

Instruction manual

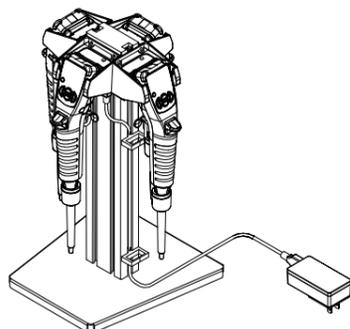

© 2018 A&D Company, Limited. All rights reserved. No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company, Limited.

The contents of this manual and the specifications of the instrument covered by this manual are subject to change for improvement without notice.



1WMPD4003079C

Safety Use

Please note the following items when the device is used.

CAUTION


- Do not disassemble the device. Any damage or loss of functions caused by disassembly is not covered by warranty.
- Do not link multiple charging stands, as it may cause fire or damage the units or the PA/PB series of electronic pipettes (hereinafter called PA/PB).
- Use the provided AC adapter, as otherwise the charging stand, the PA/PB, or the power supply devices may be damaged.

1. Introduction


This manual describes how the AX-ST-CH-M4 charging stand for four pipettes works. To fully understand and make use of the product, please read this manual carefully before use. Also, please read the instruction manual for the PA/PB.

2. Package Contents and Each Part Name

- Charging stand for four pipettes
- Stand base
- Screw (2 pcs)
- AC adapter (AX-TB283)
- Instruction manual (This document)

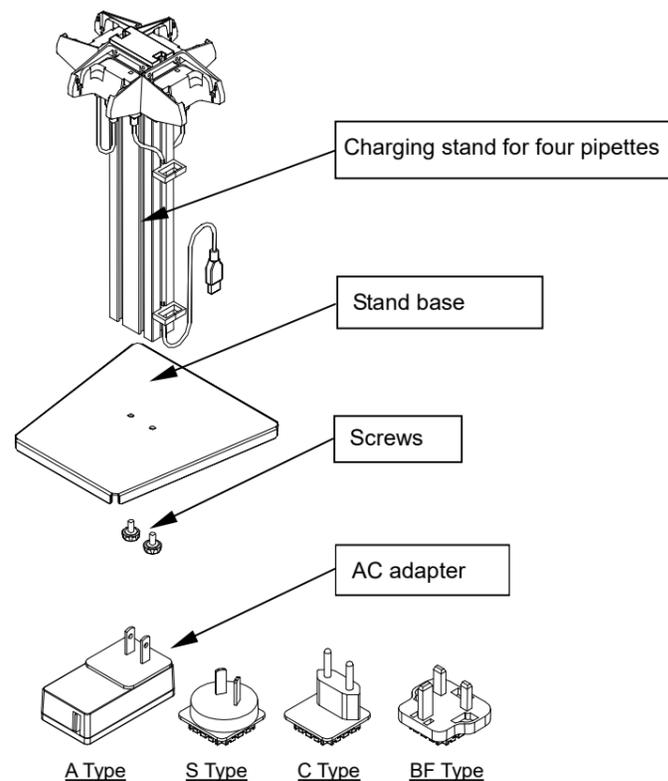

* The A-type plug is initially attached to the AC adapter for use in Japan. For use outside of Japan, please change it to the type appropriate for your country.

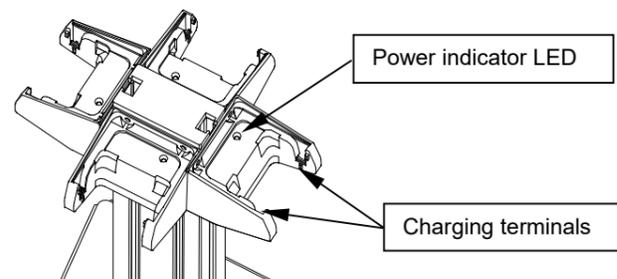

Charging terminals : Receive the PA/PB for charging.

Power indicator LED : Lights up in green when power is supplied to the charging stand.


3. Features

- Charging starts automatically when an PA/PB is hooked onto the charging stand.
- Power supply stops automatically when charging is complete.
- Space saving installation (200 mm×200 mm)
- It is possible to charge four pipettes at the same time.

4. How to Use


1. Securely mount the charging stand for four pipettes onto the stand base with the screws (2 pcs)

— Securing the post to the stand base —

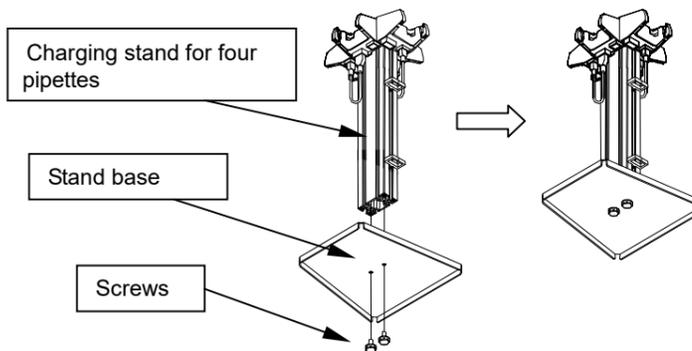


2. Prepare the provided AC adapter.

3. Connect the power cable (USB- Type A) connected to the charging stand to the AC adapter, and insert the AC adapter into an outlet. The four power indicator LEDs (Green) will light up.

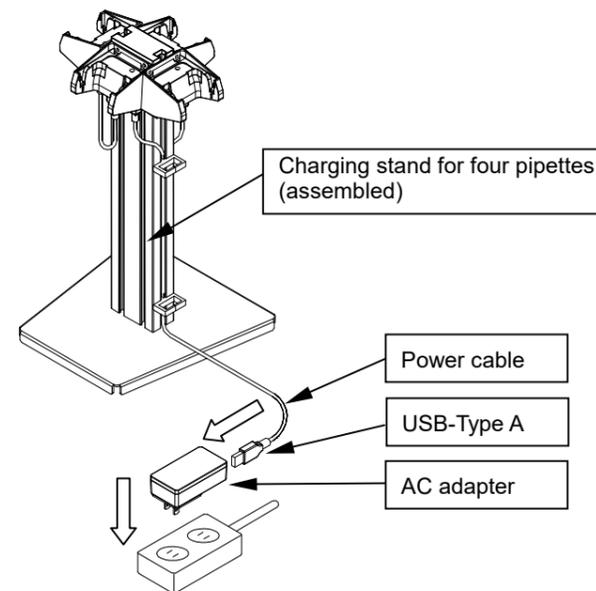

4. Hook the PA/PB onto the charging sections to start charging. Once charging starts, the battery indicator at the bottom left of the PA/PB display will blink as shown in the figure on the right.

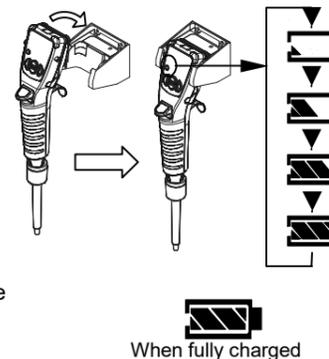

5. Charging is complete when the PA/PB battery indicator is fully lit.

6. The PA/PB can continue to be hooked to the charging stand after they are fully charged. (They will not be overcharged.)

5. Specifications

Input (Specified adapter)	100 VAC/0.3 A (50/60 Hz)
Input section / output section	DC 5V
Connectable devices	Electronic pipette PA/PB series
Dimensions	195(W)×350(H)×195(D) mm
Net weight	Approx. 1430 g (Excluding the AC adapter)

6. Accessories

- Linking cable AX-KO4730-10
- AC adapter AX-TB283