


QUICK OPERATION GUIDE

SINGLE CHANNEL ELECTRONIC PIPETTE (MPA-10 / 20 / 200 / 1200 / 10000)

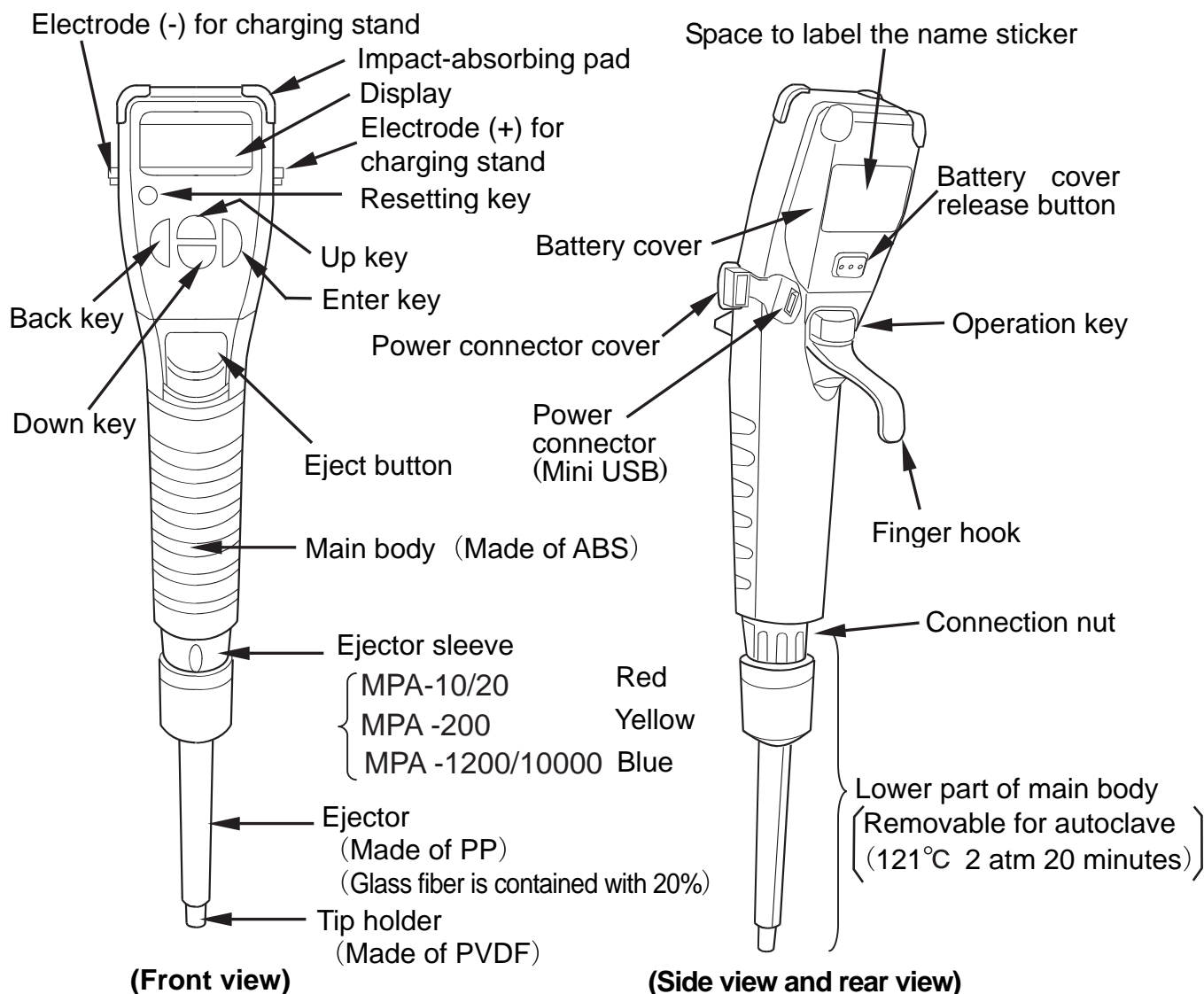


SYS System setting mode
AUTO Standard mode
MD Multiple dispensing mode
MIX Mixing mode

Aspirating and dispensing
 speed display

Battery mark

Resetting key

Back key

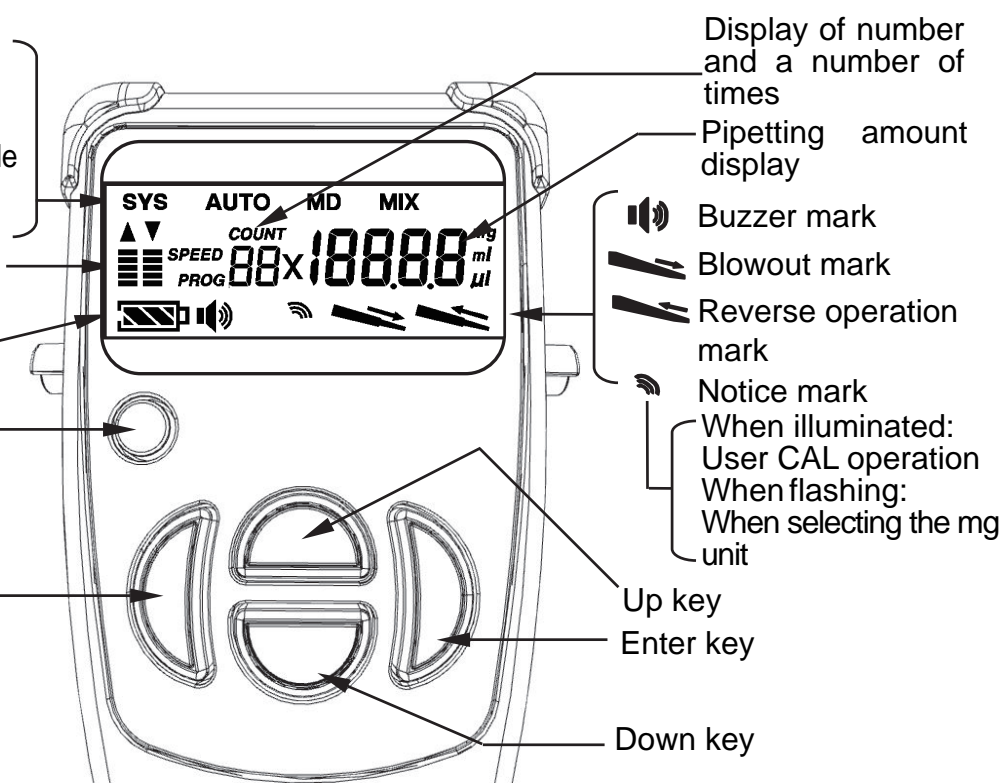


Example of the function and operation

(The following example is when using MPA-200.)


Standard mode (AUTO)

Example: When dispensing with the dispensing volume of 100µL

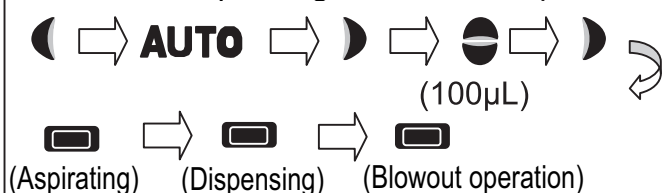


Multiple dispensing mode (MD)

Example: When multiple dispensing of 40μL
× 3 times = 120μL

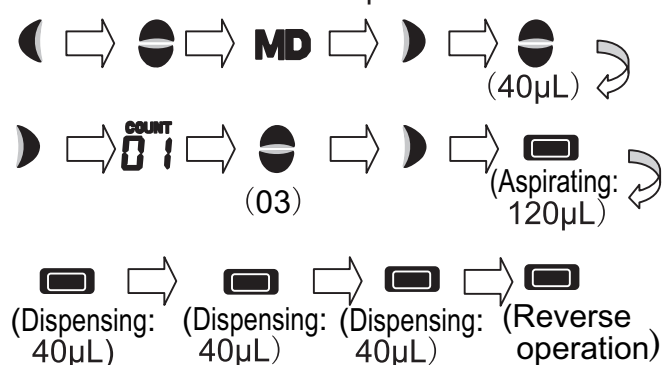


Mixing mode (MIX)

Example: When mixing with 100μL five times

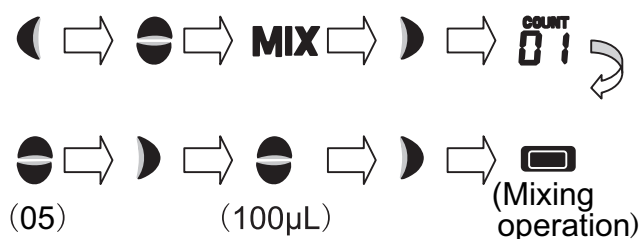


System setting mode (SYS)

Example: When setting the buzzer to OFF

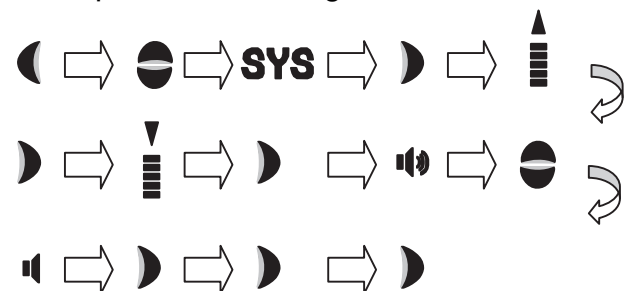


Volume calibration function (μL calibration function)

Example: When calibrating the dispensing amount to 190µL.

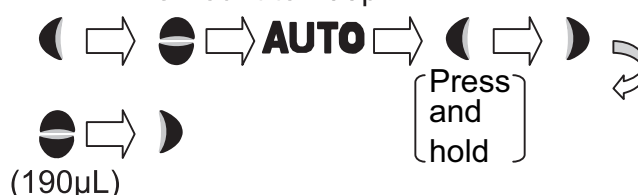

Initializing the volume calibration

Example: When initializing the dispensing amount calibrated

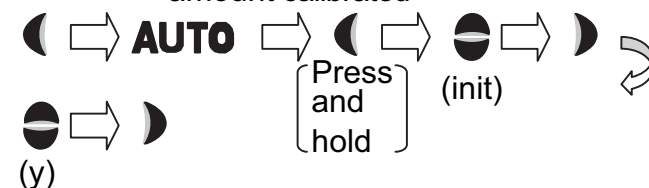


Dispensing in a unit of weight (in mg unit)

Example: when changing the unit of the dispensing amount to a unit of weight





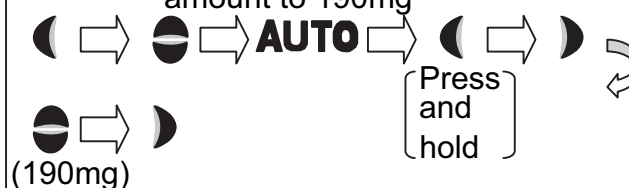

Weight calibration function (mg calibration function)

Example: When calibrating the dispensing amount to 190mg



Additional contents

In operating procedure on this page,

 means Back key,  means Up/Down key,
 means Enter key and  means Operation
 key. Press these keys when operating.