


WinCT Version Update Information

(Last updated: January 27, 2025)

2025/1/27
A&D Company, Limited.
Marketing Division

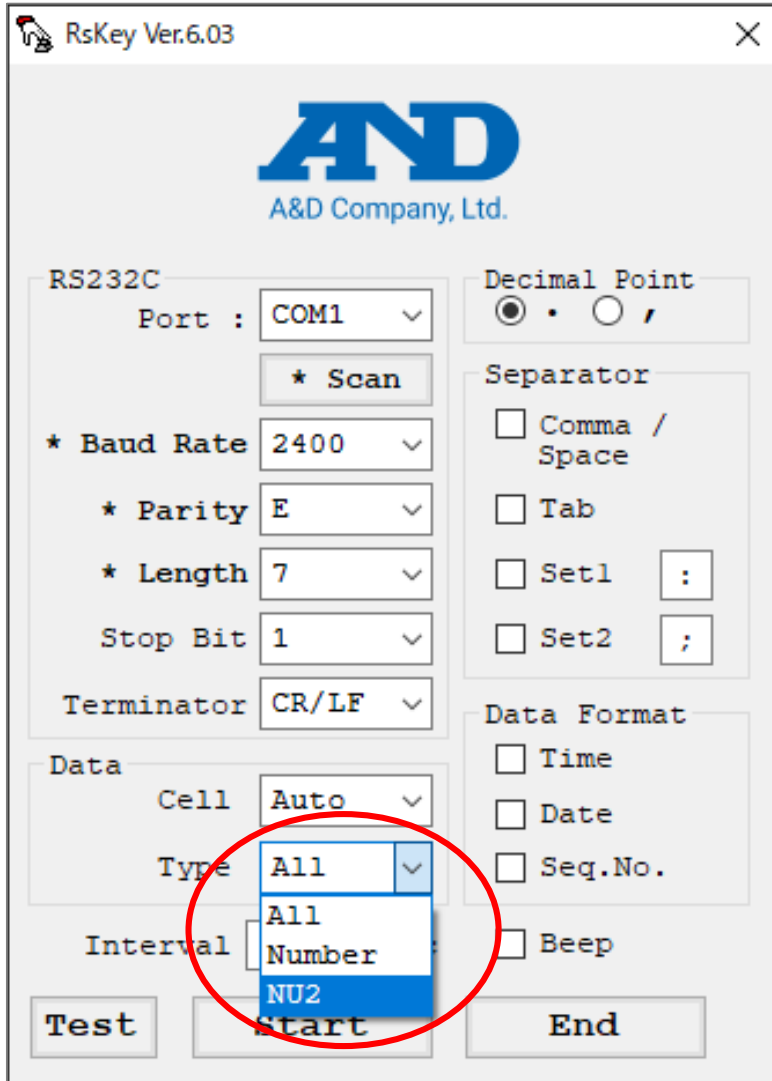

Latest Version

- Ver.6.02 >> Ver.6.03

Component Versions

- RsCom: Ver.6.02
- RsKey: Ver.6.02 >> Ver.6.03
- RsWeight: Ver.6.02

Update Details

- Added "NU2" to the output data types (Data, Type) to RsKey Ver.6,03

Details of "NU2"

- No plus sign and leading zeros (this data type is useful when the system receiving the weighing data requires the plus sign and any leading zeros to be eliminated).

Output Examples

Data Type	ALL	Number	NU2
	ST,+00162.57 g	+00162.56	162.56
	ST,-00002.57 g	-00002.57	-2.57